



Uppsala December 4th, 2025

Release Notes

Olink Target Panel Data Archive for NPX™ Signature software v2.0 onwards

Release version: OlinkTarget_PanelDataArchive_2025NOV27_NPXsv2.dat

Release Date: 2025DEC04

Included in Olink Target Panel Data Archive 2025NOV27

Added Panel	Content
Olink® Target 48 and Olink® Target 96	Support for ALL previously released Olink® Target 48 and Olink® Target 96 reagent batches.
Newly added Olink® Target 48 Panels	A new Panel Data File for Olink® Target 48 Cytokine is available, Panel Data File version: 932012.
Newly added Olink® Target 48 Panels	A new Panel Data File for Olink® Target 48 Mouse Cytokine is available, Panel Data File version: 934101.

Note Olink® Target 48 Cytokine:

In Target 48 Cytokine kits connected to panel data file version 932012 a new antibody is included for the IL17F assay. As this is a new antibody the assay has gotten a new OID to highlight this.

Note Olink® Target 48 Mouse Cytokine:

The panel data file for the new Target 48 Mouse Cytokine reagents, panel data file version: 934101 contains updates that are not included in previous panel data file versions (panel data files 934001 & 934002). These panel data file updates are not compatible with the two previous reagent lots.

www.olink.com

© 2025 Olink Proteomics AB, part of Thermo Fisher Scientific.

Olink products and services are For Research Use Only. Not for use in diagnostic procedures.

This document is not intended to convey any warranties or representations of any kind.




In NPX Signature v2.0.2 it is not possible to change the panel version between the new (934101) and the two previous ones (934001 & 934002) in an already started project, as this will lead to an application crash. If the wrong panel data file version is mistakenly chosen it is needed to create a new project.

In NPX Signature v2.2 it is possible to change the panel version between the new (934101) and the two previous ones (934001 & 934002) in the project, if the project is created in NPX Signature v2.2.

In NPX Signature v2.1 and NPX Signature v2.2 (for projects created in previous software versions) it is not possible to change the panel version between the new (934101) and the two previous ones (934001 & 934002) in the project, an error message will occur. If the wrong panel data file version is mistakenly chosen it is needed to create a new project.

Note Manage Panel Data Files:

In NPX Signature v2.2 there is a known issue that after uploading the new panel data archive the name of the newly uploaded and previously uploaded panel data archive is displayed at the same time:

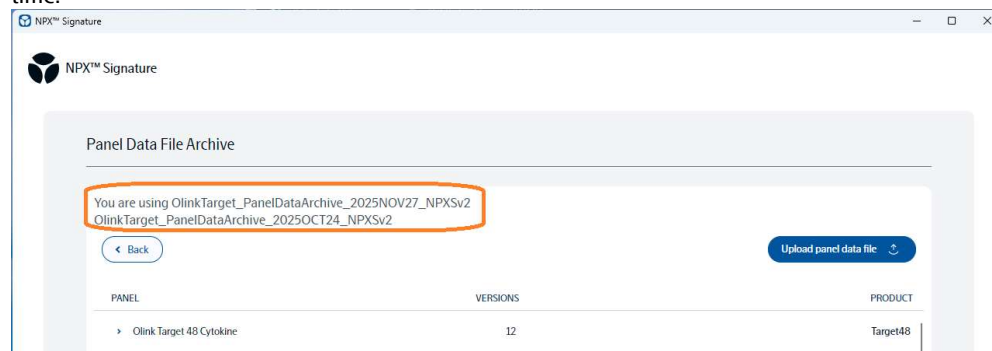

This is a visual issue only, panels and panel version are displayed and can be used correctly.

Software download

The latest version of NPX Signature software and the Target Panel Data Archive is available to download from <https://olink.com/software/download>

Password: "bg7Y0??yg89&!"